

**Rely+On™
Virkon™**

An important message from the manufacturer:

Rely+On™ Virkon™ highly effective against Human Corona Virus

**Inactivation of the Human Corona Virus:
Rely+On™ Virkon™**

A recent study shows that Rely+On™ Virkon™ is highly effective at inactivating the Human Corona Virus.

Please read more on the following message from Anneliese Bischof, Business Director Disinfection, Lanxess



Rely on Virkon is distributed in New Zealand by Bio-Strategy Ltd

Bio-Strategy also supply wipes that when impregnated with a 1% solution of Rely+On™ Virkon™, offer a convenient and consistent way to apply this high level disinfection to hard surfaces.


Preparation of wipes

Safety notice: Gloves and goggles must be worn.

- A. Prepare a 1% Virkon solution. Fill a suitable container with 500ml warm water and add 1 x 5g Virkon Tablet. Allow to dissolve fully.
- B. Open wipe pack RHS V1, carefully pour (reconstituted) Virkon solution into the pack opening.
- C. Seal pack and rotate until all wipes are saturated with VIRKON liquid.
- D. In space provided, write on the date the mixture was prepared.
- E. Use wipes as directed until container is empty or expiry date reached.



**minimize the risk for yourself and/or the team
call 0800 34 24 66 to order**

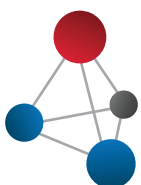

bio-strategy
delivering technology

VIRKON, SUPPLIED BY:
Bio-Strategy Limited | T: 0800 34 24 66
sales.nz@bio-strategy.com | www.bio-strategy.com

Rely+On™ Virkon is the registered trade mark of **LANXESS**

Antec International Limited, Sudbury, Suffolk CO10 2XD, United Kingdom

24 January 2020

Rely⁺On™ Virkon™ highly effective against Human Coronavirus

As you are aware, Human Coronavirus continues its rapid spread across Europe, China and Southeast Asia, with suspected cases in the UK and Canada. Many authorities have announced screening measures for passengers from China, including the major airport hubs of Dubai and Abu Dhabi.

With no vaccine or other medical treatment currently approved, we are pleased to advise that Rely⁺On™ Virkon™ has already received independent evaluation against the virus.

Rely⁺On™ Virkon™ met the performance requirements specified in the study protocol. The results indicate complete inactivation of Human Coronavirus under these test conditions, as required by the U.S. EPA. Tests were carried and confirmed by independent company ATS labs in 2015.

Rely⁺On™ Virkon™ achieved complete inactivation of the Human Coronavirus at a 1:100 dilution rate with a 10 minute contact time.

These test results provide independently proven data on the ability of Rely⁺On™ Virkon™ to effectively and completely kill the Human Coronavirus rapidly.

Yours sincerely



Anneliese Bischof
Business Director Disinfection

Antec International Ltd
Material Protection Products
Windham Road
Chilton Industrial Estate
Sudbury, Suffolk CO10 2XD
United Kingdom
01787 377305

www.lanxess.com
www.virkon.com

Board of Management:
Jacques Louis Albert Perez
Christopher James Flett
Dr. Juan González-Blanco

Registered Office:
Windham Road
Chilton Industrial Estate
Sudbury, Suffolk CO10 2XD
No. 690279 England